


TBSI 5-CH Headstage to FHC electrode Adapter Instructions for Use

PN 7118

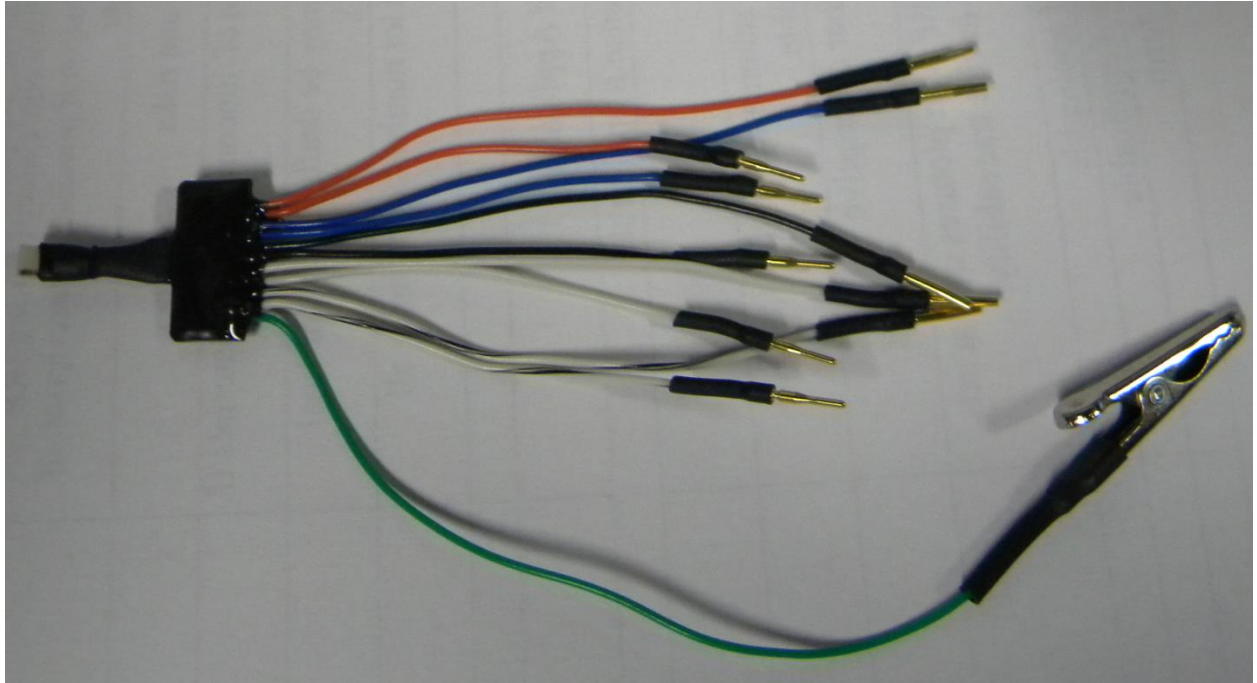

Blackrock Microsystems, Inc.

391 Chipeta Way, Suite G

Salt Lake City, UT 84108


Tel: (866) 806-3692

www.blackrockmicro.com


TBSI 5-CH headstage to FHC electrode Adapter Instruction for Use

1 Hardware

1.1 Blackrock TBSI 5-CH headstage to FHC electrode Adapter Pin-out

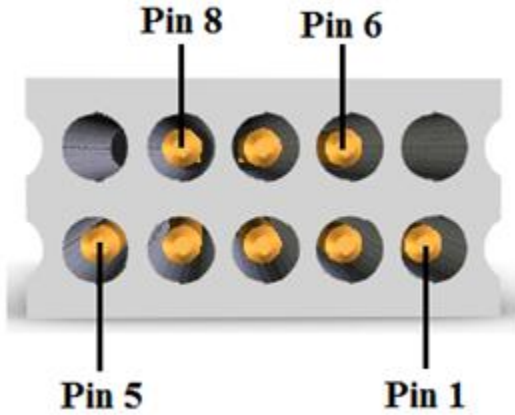

Pin	Pin Connection
1	Channel 5
2	Channel 4
3	Channel 3
4	Channel 2
5	Channel 1
6	VSS
7	ACGND (Animal Ground) / REF

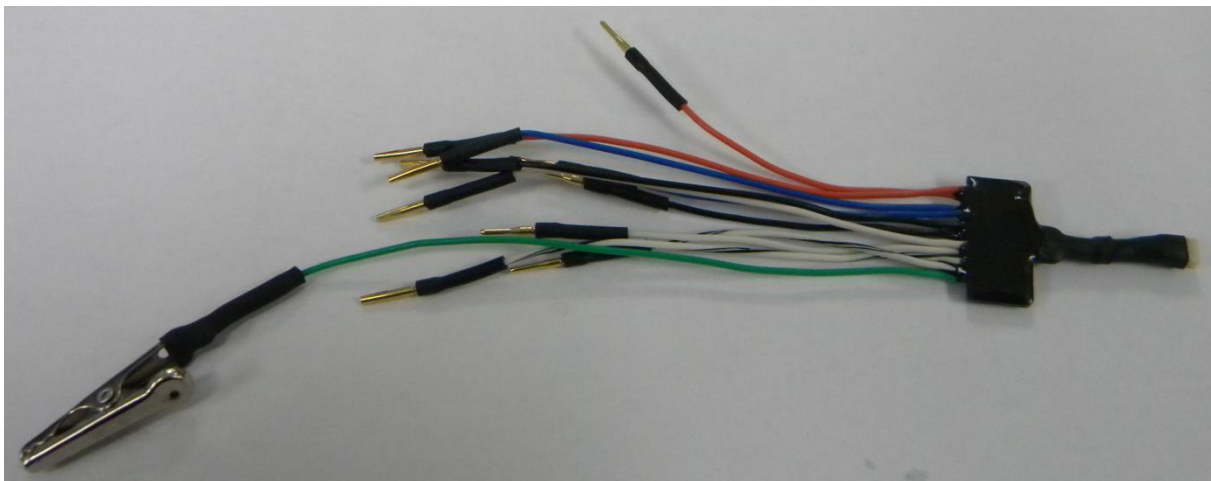

Male Pins (GND)			Female Pins (Electrodes)		
Wire Color	Pin #	Connection	Wire Color	Pin #	Connection
Orange	2	GND	Orange	1	Electrode 1
Blue	6	GND	Blue	5	Electrode 2
Black	10	GND	Black	9	Electrode 3
White	14	GND	White	13	Electrode 4
White w/Black	18	GND	White w/Black	17	Electrode 5
Green w/Alligator Clip	22	ACGND (Animal GND) / REF			

Figure 1-2: TBSI 5-CH headstage to FHC electrode Adapter Pin-Out

TBSI 5-CH headstage to FHC electrode Adapter

Instruction for Use

1.2 Adapter Ground Connections

The Omnetics FHC-Style Adapter is able to handle up to 5 electrodes at one time. Any un-used Electrode connection should be shorted to ground by mating it with the corresponding ground pin.

2 Warranty

Blackrock Microsystems, Inc. warrants that its products are free from defects in materials and manufacturing for a period of one year from the date of shipment. Blackrock will, at its option, repair or replace any product that does not comply with this warranty. This warranty is voided by:

1. Any modification or attempted modification to the product done by anyone other than an authorized Blackrock employee
2. Any abuse, negligent handling or misapplication of the product.

This constitutes the sole warranty made by Blackrock, Inc. There are no other warranties, expressed or implied, which extend beyond those described herein or to anyone other than the original purchaser, including the implied warranties of merchantability and fitness for a particular purpose. In no event shall Blackrock Microsystems, Inc. be liable for any incidental or consequential damages, or for the infringement of any patent rights or third party rights due to the use of its products.

2.1 Return Merchandise Authorization (RMA)

In the unlikely event that your Adapter needs to be returned to Blackrock for repair or maintenance, do not send any equipment back without a Return Merchandise Authorization Number. An RMA number will be issued to you by a Blackrock representative. If you need to obtain an RMA number, you may contact a product support representative at (801) 582-5533 or toll free at (866) 806-3692.

Once an RMA number has been issued, it is important to safely pack the returned item for shipping back to Blackrock. It is preferred that you save the original boxes and packing materials that your system arrived in for return shipment. Please address the package as follows:

Blackrock Microsystems, Inc.
ATTN: RMA#
391 Chipeta Way, Suite G
Salt Lake City, UT 84108 USA
Tel: (801) 582-5533