


Custom Commutator to HDMI Adapters

PN-8849 & PN-8850

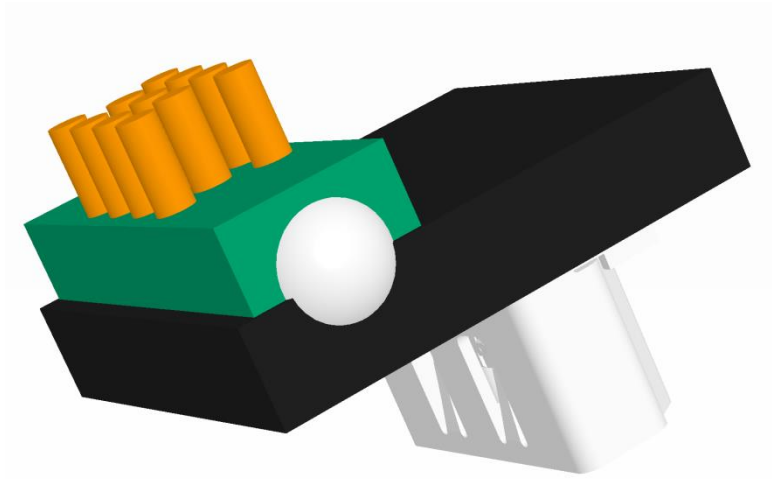

Blackrock Microsystems, LLC

630 Komas Drive • Suite 200
Salt Lake City • UT 84108

T: +1 (801) 582-5533


www.blackrockmicro.com
support@blackrockmicro.com

1 Introduction


The adapter provides interface between HDMI cable for Cereplex headstages and a commutator with 7-Pin Dale connector and DB-15 connector.

2 Hardware

2.1 Blackrock 7-Pin Dale – HDMI-A Pin-out PN-8849

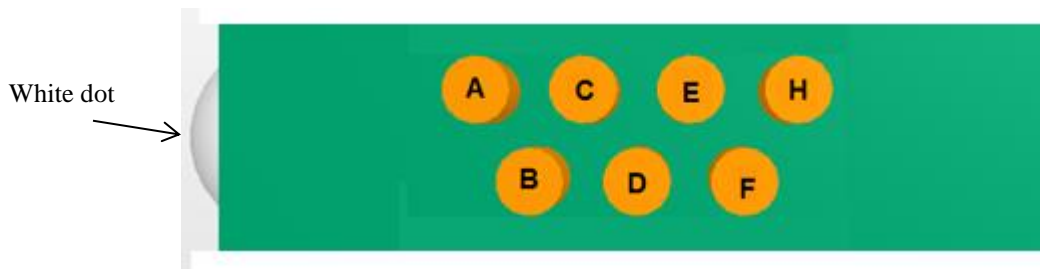


Dale MMP22G5-7

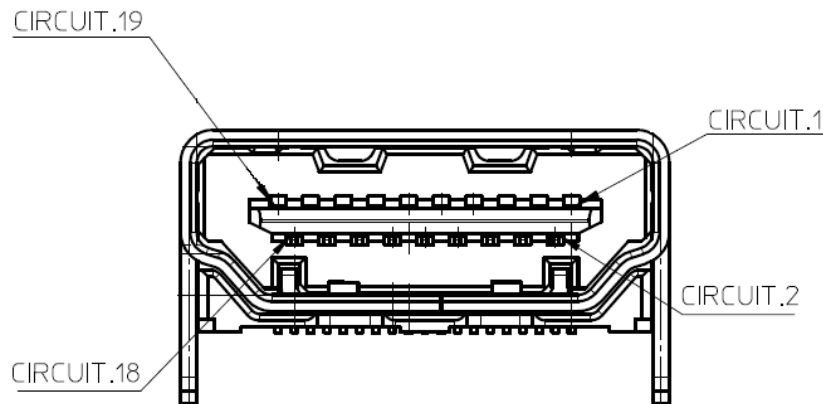


HDMI Molex 04659-1000

DALE	SIGNAL	HDMI
C	GND	2, 5, 8, 11, 19
F	+V	18
H	-V	17
A	+CLK	10
D	+DATA	7
E	-DATA	9
B	-CLK	12

2.2 Blackrock DSUB 15-Pin – HDMI-A Pin-out PN-8850

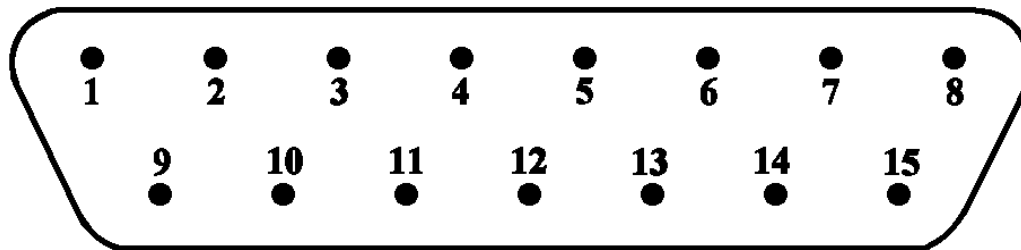

DSUB-15 5-747908-2

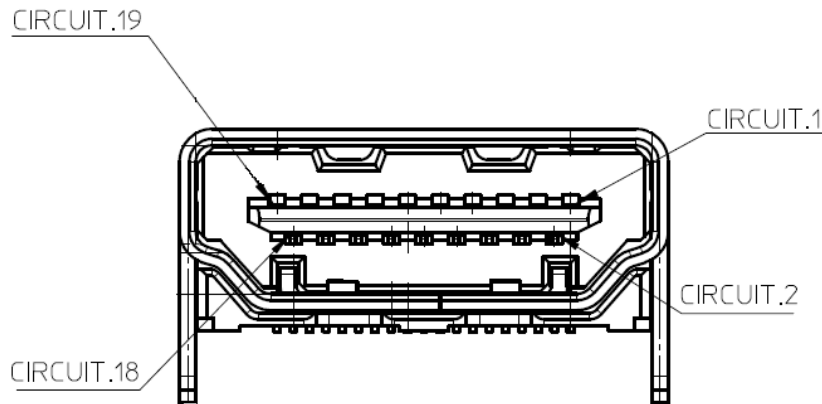

HDMI Molex 04659-1000

DALE	SIGNAL	HDMI
3	GND	2, 5, 8, 11, 19
6	+V	18
7	-V	17
1	+CLK	10
4	+DATA	7
5	-DATA	9
2	-CLK	12

3 Warning

These assemblies can only be used for animal research/applications.

4 Warranty

Blackrock Microsystems warrants that its products are free from defects in materials and manufacturing for a period of one year from the date of shipment. Blackrock will, at its option, repair or replace any product that does not comply with this warranty. This warranty is voided by:

1. Any modification or attempted modification to the product done by anyone other than an authorized Blackrock employee
2. Any abuse, negligent handling or misapplication of the product.

This constitutes the sole warranty made by Blackrock, There are no other warranties, expressed or implied, which extend beyond those described herein or to anyone other than the original purchaser, including the implied warranties of merchantability and fitness for a particular purpose. In no event shall Blackrock Microsystems, be liable for any incidental or consequential damages, or for the infringement of any patent rights or third party rights due to the use of its products.

4.1 Return Merchandise Authorization (RMA)

In the unlikely event that your adapter needs to be returned to Blackrock for repair or maintenance, do not send any equipment back without a Return Merchandise Authorization Number. An RMA number will be issued to you by a Blackrock representative. If you need to obtain an RMA number, you may contact a product support representative at (801) 582-5533 or toll free at (866) 806-3692.

Once an RMA number has been issued, it is important to safely pack the returned item for shipping back to Blackrock. It is preferred that you save the original boxes and packing materials that your system arrived in for return shipment. Please address the package as follows:

Blackrock Microsystems
ATTN: RMA#
630 Komas Drive, Suite 200
Salt Lake City, UT 84108 USA
Tel: +1 (801) 582-5533